Transmission Electron Microscopy

F90 - Fortgeschrittenen Praktikum UNIVERSITÄT HEIDELBERG

> Anne Kast, Rasmus Schröder Revised: Wolfgang Köntges

> > September 20, 2019

Important Information:

Please carefully read the email you receive upon booking this experiment for further information. We recommend to use the literature given in the link for course preparation. The literature contains the neccessary theory of operating a transmission electron microscope.

Section 3 will help you in your prepartion especially to find the necessary information from the literature. The tasks intend you to think about and discuss the main topics of this course. Perfect answers and correct solutions in every detail is not necessary to attend this course successfully.

Furthermore we recommend to use the image program Fiji for your analysis. Notice the information in appendix B concerning Fiji.

Contents

1 I	ntroduction	1
2 C	Outline of Experiments	5
2.1	Day 1: Introduction, Basic Operation and First Experiments	5
2.2	Day 2: Diffraction, Bright- and Darkfield Imaging	5
2.3	Day 3: Elective Experiments TEM or SEM	6
3 C	Course Preparation	7
3.1	The Imaging system of a TEM	7
3.2	Contrast Transfer Function	8
3.3	Diffraction, Bright- and Darkfield-Imaging	9
3.4	Scanning Electron Microscopy (SEM)	9
3 E	Expectation of Course Preparation	13
3.1	The Imaging system of a TEM	13
3.2	Contrast Transfer Function	18
3.3	Diffraction, Bright- and Darkfield-Imaging	22
3.4	Scanning Electron Microscopy (SEM)	23
4 E	Experimental Procedure	25
4.1	Day 1: Getting Started	25
4.2	Day 2: Diffraction Experiments	30
4.3	Day 3: Elective Experiments at the TEM or SEM	35

ii	Contents							
4.3	B.1 Elective Experiments at the TEM	35						
4.3	B.2 Elective Experiments at the SEM	38						
5 E	xpected Documentation of the Experiments	39						
5.1	Magnification Calibration	39						
5.2	Judging Image Quality	40						
5.3	Diffraction, Bright- and Darkfield Imaging	40						
5.4 Understanding the Contrast Transfer Function								
5.5	Investigating Vegard's Law	41						
5.6	6 Comparison between TEM and SEM							
Bibli	ography	43						
Appe	endices	45						
A D	iffraction Gold	47						
вн	elpful Features in Fiji	49						
B.1	Opening images and scalebars	49						
B.2	Adjusting Brightness and Contrast	49						
B.3	Fast Fourier Transforms (Power Spectrum)	50						
B.4	Image Stacks	50						
B.5	Radial and Line Profile	50						
C P	rovided Samples	53						

1. Introduction

Transmission electron microscopy has in recent years developed into a – somewhat – versatile tool for physical as well as biomedical sciences. The basic interaction between electrons and sample, i.e. the interaction with the atoms in the sample consisting of atomic nuclei and electrons, is well understood and can be described by theories for the elastic electron-nucleus, and the inelastic electron-electron interaction (cf suggested literature, e.g. Williams and Carter [Wil09], Reimer and Kohl [Rei08], R. Erni [Ern10], J. Rodenburg [Rod18]). However, necessary instrumentation for high resolution, atomic imaging in a transmission electron microscope (TEM) had to be developed in a long process culminating in the correction of lens aberrations realized in the last two decades. The correction of lens aberrations such as spherical aberration, coma, and more recently also chromatic aberration has allowed scientists to directly visualize columns of atoms in crystalline material and also individual carbon atoms in graphene [Hai98]. Figures 1.1 and 1.2 were adopted from the SALVE web-page (http://www.salveproject.de), a recent collaborative project for optimizing imaging and electron spectroscopy on carbon-based materials.

In material science today, the TEM is used for high-resolution imaging as illustrated here, but also for studies based on more conventional electron scattering experiments, i.e. electron diffraction on crystalline material, as well as for analytical studies. Such studies utilize the special nature of inelastic interactions between the incident electrons and the electrons in the sample. In general, all these techniques are employed to obtain information about the structure of a material and its localized electronic



Figure 1.1: Strontium titanate as a typical materials science sample where in the past it was not possible to image individual columns in an oriented and thinned specimen. (A) shows the known atomic structure, which has been imaged in the past in conventional microscopes as shown in (B). Such images could only be interpreted by advanced simulation of image contrast from known models. (C) and (D) show images and more modern simulations of aberration corrected imaging. Figure adopted from SALVE web-page.



Figure 1.2: Examples of conventional (upper row) and aberration corrected images (lower row) of a mixed crystalline/amorphous Si specimen. The direct imaging in the aberration corrected microscope shows the individual Si atom columns of the crystalline area (left part of the picture). This cannot be resolved in the conventional image. Figure adopted from SALVE web-page.

properties. Imaging resolution is very often in the 1 Å to even sub-Å range. The resolution of the analytical signal is normally reduced to the typical delocalization of the inelastic interaction in the sample.

In biomedical electron microscopy, the situation is quite similar: With advances in sample preparation and, in particular, constantly improving electron imaging detectors, it is now possible to image biomolecules such as e.g. proteins or protein complexes in their 3D structure at a quasi molecular level, sometimes in the range of 3 to 5 Å – more routinely in the range of 6 to 8 Å. Figure 1.3 shows a typical image of such a «6 Å study» of the actin-myosin model complex, i.e. the force producing protein complex, which works in our muscles.

If one compares the images of a materials sample with the images of the actin myosin complex one clearly realizes, that the level of resolution and atomic detail is completely different. The reason for this is very simple: The electrons used for imaging also damage the sample. This is obvious from the physics of inelastic interactions, their energy transfer and the fact, that inelastic interactions are more probable than elastic interactions for light material. For large Z materials (Z > 26), elastic interactions start to win. Biomedical samples, which are prepared in a glassy, vitried



Figure 1.3: (A) Images acquired by transmission electron cryo microscopy of F-actin decorated with myosin Va motor domains with truncated lever arm, in nucleotide-free Rigor state. The defocus value is $1.7 \,\mu m$ and $5 \,\mu m$ for the insert, respectively. (B) The reconstructed density map of the nucleotide-free complex shows the typical helical symmetry of the filamentous actin-backbone. Shown are six actin molecules in their ribbon representation (light yellow) and four myosin motor domains (blue). Images from our own paper [Wul16].

aqueous buffer layer of about 50 to 100 nm thickness, are simply «boiled away» by our electron beams. Thus, images have to be taken at lower magnication and lower electron dose, which decreases resolution and increases noise.

What will you learn from this experiment?

As you might expect, TEMs which deliver sub-Å resolution or allow the imaging of a frozen sample at liquid nitrogen temperature are very specialized, expensive, and most of the time used for our research. A microscope – as small as the one used here in our practical course – will not have all this expensive instrumentation, but will get you acquainted with some very important physical facts, which we in our daily research still have to be aware of. Familiar topics in this respect are the optimal optical alignment of an imaging system in particle optics, basic imaging properties when working with a particle beam, the recording of an image and basic image processing steps. Also, electron diffraction experiments can be performed with our microscope, and very fundamental experiments to test the coherence of an electron beam and how it can be inuenced by different beam forming parameters. In the end you should have a first feeling – and also the physical basis for it – what it takes to set up a microscope optimally and how to tune it for highest performance. One of the «holy grails» of transmission electron microscopy, the Contrast-Transfer-Function (CTF), should then be something, which has lost its legendary aura. While we provide a guideline for experiments with our samples (which we provide as well), we also will give you samples for «free microscopy». This will be tissue samples from biological specimens, plant root or mouse muscle prepared with different protocols. Besides all the proposed experiments, documentation, suggested reading, and discussion, we would like to encourage you to simply «play around» with these samples – following the rule that «one can see a lot just by looking». To make this worthwhile – however – we will discuss your images (and your documentation of this free microscopy) with you and also explain a bit of the biology of the samples.

2. Outline of Experiments

2.1 Day 1: Introduction, Basic Operation and First Experiments

Theory and basic operation of the transmission electron microscope (TEM)

- Experiment 1.1: Acquire a defocus series of the myelin sheath in a mouse muscle. First insights into the effect of the Contrast Transfer Function.
- Get to know the lens system of the microscope.
- Experiment 1.2: Determine the effect of a) over- and underfocus, b) astigmatism, c) coherence, d) centered objective aperture and e) different magnification settings on the image quality.
- Removement and insertion of the sample holder. Alignment of the microscope.
- **Magnification Calibration**: Record images of a standard sample for magnification calibration.

2.2 Day 2: Diffraction, Bright- and Darkfield Imaging

Diffraction Experiments and Basic Principles of Bright- and Dark-Field Microscopy

- Understand the formation of diffraction patterns and which information they provide about the sample.
- Experiment 2.1: Acquire a diffraction pattern of a thin film (Au) at 150 mm or 250 mm camera length.
- Understand different imaging methods in the TEM using the diffraction mode.
- Experiment 2.2: Acquire bright- and darkfield images of MgO crystals.

Elective Experiments for TEM or SEM

You can choose one of the following experiments depending on what you want to do the next day:

- Experiment 2.3: Record a second DP of the gold film at another camera length for calibration purposes.
- Experiment 2.4: Take images of one or two provided biological samples to compare the images with the according SEM images.

2.3 Day 3: Elective Experiments TEM or SEM

You can either continue with experiments at the TEM or the SEM.

Further Experiments at the TEM (diffraction, bright- and darkfield imaging

- Experiment 3.1: Vegards law: Acquire diffraction patterns of alloys (Au/Pd or Pt/Pd) at the camera lengths you used the previous day.
- Experiment 3.2: Acquire a defocus series of a catalase crystal to further understand the effect the Contrast Transfer Function on the image.
- Experiment 3.3: Measure and use the periodicity of a catalase crystal to calibrate higher magnifications of the TEM.

Elective Experiments at the SEM (biological samples)

- Get to know the basic operation of a SEM (magnification, focus, wobbling, astigmatism correction).
- Experiment 3.4: Acquire images of samples that you already investigated with the TEM for comparison.
- Understanding the differences between TEM and SEM. Explaining advantages/disadvantages of both methods.
- Experiment 3.5: Acquire images of biological samples you can't image in the TEM.

3. Course Preparation

Work on the following tasks and be prepared to discuss them over the course of the three days.

3.1 The Imaging system of a TEM

The beam paths in electron optics can be drawn analogous to light optics. Chapters 6 and 9 of [Wil09] may help you working on the following tasks.

- 1. Name reasons, why we are using microscopes.
- 2. Consider **two** scattering points and **one** non-scattering point (vacuum) located in front of an ideal lens as depicted in figure 3.1 at the end of this section. You can use the figure for the following constructions.
 - Define the image plane of the lens at an arbitrary, but reasonable place behind the lens (arbitrary magnification).
 - Construct the image of the given objects by using the centre of the lens and the previous defined image plane. You may benefit from your knowledge of optical geometry.
 - Assume that all electrons (scattered and non-scattered) passing a point are moving through the according image point in the image plane as well. With this construct the beam paths of non-scattered (0°) and scattered (α) electrons at the following angles α in different colors: 40°, 20°, -20°, -40°
 - Mark characteristic points and planes on basis of your constructions and explain, why they are characteristic.
 - Mark in the same figure the location of the following apertures and explain their purpose.
 - Condenser aperture (CA)
 - Objective aperture (OA)
 - Selected-area aperture (SA)
- 3. Explain the purpose of the following parts in a basic TEM setup: Condenser, Objective, Projective.

3.2 Contrast Transfer Function

Sections 6.5 and 28.7 of [Wil09] may help you working on the following tasks.

- 1. Explain «contrast».
- 2. Explain the difference between the following contrast mechanisms and indicate the ones used for imaging in a conventional TEM.
 - absorption-contrast
 - aperture-contrast
 - phase-contrast
 - filter-contrast
- 3. Explain in a few sentences the difference between «Fresnel» and «Fraunhofer» diffraction.
- 4. Name possible lens aberrations and how they can be compensated.
- 5. Explain «Scherzer-Focus» and its significance. (No calculation needed! This can be found in section 28.7 of [Wil09])
- 6. You will have to calibrate all magnifications of the microscope in real space and fourier space of an image showing a standard sample. One calibration method is described in Chapter 9 of [Wil09] (p. 164f).
 - Download the image «Calibration.tif» from the link showing an image of a standard specimen so called «cross grating» (2160 lines/mm) with latex beads on it (black dots with \emptyset 0.261 µm).
 - Open the image in an image program (e.g. Fiji) and get an idea, how to measure lengths and intensity profiles with this program. You will need that for data analysis later on. Helpful features of Fiji can be found in appendix B.
 - Find out what measurements need to be done in real space **and** in the fourier transformation of the image (distance real, diameter latex beads, distance FFT) to calculate the pixel size in the image.
 - Argue, why the magnification is calibrated, if you know the pixel size of all magnification settings and estimate the error you would expect from your measurements.

3.3 Diffraction, Bright- and Darkfield-Imaging

Chapters 11 and 12 of [Wil09] may help you working on the following tasks.

- 1. Write down the «Bragg condition».
- 2. Describe, what kind of diffraction pattern you would expect from a single crystal and therefore for a polycrystalline sample.
- 3. Write down «Vegard's law» and calculate the lattice parameter for Pt/Pd and Au/Pd (each 80/20).
- 4. You will have to calibrate the camera length of the microscope. As for the previous calibration a suitable method is described in Chapter 9 of [Wil09] (p. 165f).
 - Download the image «DiffCalib.tif» from the link. This is an image stack containing 8 images of a polycrystalline gold diffraction pattern.
 - Open the image stack in an image program (e.g. Fiji) and use the «Z-Project» function of Fiji (see appendix B) since this will help you to present your data later on.
 - Use the «Radial Profile» function in Fiji to calibrate the camera length. Estimate the error, you would expect from your measurements.

3.4 Scanning Electron Microscopy (SEM)

The following preparation part only needs be done, if you choose the elective experiments for the SEM. You can do after the second part of the course, after you decided to go for the SEM experiments. Chapters 1 of [Rei93] may help you working on the following tasks.

- 1. Name the principle differences between TEM and SEM.
- 2. Explain the image formation in the SEM in contrast to the TEM.
- 3. Describe qualitatively the interactions between the primary electron beam and the sample, that are essential for imaging in the SEM.
- 4. Name the resulting electrons from these interactions, which can be detected in the SEM.
- 5. Name the limiting factors for the image resolution in the SEM and compare it with the resolution limit in the TEM.

3. Course Preparation



ideal lens

optical axis

Figure 3.1: Construction of scattered and unscattered beam paths through an ideal lens. The filled dots in front of the lens represent scattering objects, while the small circle represents a non-scattering object (vacuum). The objects are illuminated homogeneusly by coherent electrons, that move parallel to the optical axis of the lens. The cross on the lens marks its center.

3. Course Preparation

4. Experimental Procedure

Always make notes of the imaging conditions at which you acquire data/images. Since you want comparable data or be able to reproduce it in the future, this is a very important experimental practice. For this purpose a spreadsheet template is provided, where you can enter all important imaging parameters. Decide who operates the microscope and who enters the microscope parameters. You can switch during the course.

4.1 Day 1: Getting Started

All important components (apertures, valve V3,...) and knobs are indicated in the figures 4.1 and 4.2.

- 1. The microscope will be turned on and adjusted by the instructor for the first experiment.
- 2. Let the instructor show you the basic operation of the microscope. During this course you will understand, what the single knobs are doing in the microscope.

The aim of the first experiments is to see the influence of the microscope settings on the image quality. We will start with the focus of the objective lens.

Experiment 1.1: Visualizing the myelin sheath of a mouse muscle

- 1. Acquire a so called defocus series of the myelin sheath in a mouse muscle and take a close look at the membranes of the myelin. Use a focus range given by the instructor in steps of 0,1 and **describe** significant differences in the images.
- 2. Compare the power spectrum (FFT) of all taken images and describe, how it changes. Do yo recognize the «Thon Rings»? Discuss these with the instructor, what the power spectrum already tells you about the image quality. You should be able to explain the influence of the focus settings on the image in real and fourier space.
- 3. For your documentation: Save the images as stack. Don't forget to note the microscope parameters in an excel sheet. Save the template sheet as a new file containing the actual date and your name.



Figure 4.1: The ZEISS EM10EL. Condensor aperture (CA), objective aperture (OA), selected-area aperture (SA), as well as important components are indicated.

Experiment 1.2: Influence of different microscope settings on the image quality

- a) The effect of over- and underfocus
- 1. The projective of the microscope is set at an magnification around 20 kx. Make sure the camera is removed from the beam path and change to a smaller magnification of 400 x (Low MAG).
- 2. Look at the fluorescent screen in low-mag-mode. You should be able to identify the embedded muscle sample.
- 3. Go to a sample area without the resin sample, in which the muscle is embedded, containing only holes of the carbon holey film .
- 4. Go back to a higher magnification of 30 kx, activate the «rabbit mode» of the camera and carefully insert the camera. You should now see an image of the holey carbon film.
- 5. Center a small hole (it should fit well on the camera) with the goniometer using the camera in «rabbit mode».

- 6. Toggle the **FOCUS** knob and observe the image.
- 7. Describe how to determine from the image (Fresnel Fringes) and the power spectrum (Thone Rongs) that your sample is in focus and adjust the microscope accordingly.
- 8. For your documentation: Make sure you have around 6000 counts and take images of holes in focus, over- and underfocus. Explain how you can discern them.

b) The effect of astigmatism

- 1. Make sure your sample is in **focus** as you learned in the previous step.
- 2. Set one of the objective stigmators (**OBJ STIG**) to maximum, which induces astigmatism to the image. **Describe how the image changed.**
- 3. For your documentation: Record an image (6000 counts) of a hole, which shows astigmatic characteristics. Explain the effect by describing the different appearance of the «Fresnel» fringes in the image.
- 4. Reset the Objective Stigmator (**OBJ STIG RECALL**).

c) The effect of coherence

- 1. Make sure your sample is in **underfocus**, so that you can see the «Fresnel» fringes.
- 2. Record an image using a small **SPOTSIZE** and short exposure times. Be careful **not to overexpose** the camera but make sure you have around 6000 counts. If neccessary increase the magnification.
- 3. Increase the **SPOTSIZE** (values of 20.0 or higher) and take an image. You will need a longer exposure time, since the countrate will drop. You might not reach the 6000 counts, not even at 7500 ms.
- 4. For your documentation: Describe the differences between images at small and high SPOTSIZE. Explain them and take the «Fresnel» diffraction into account.

d) Aperture contrast

1. Increase the sample illumination (go back to a smaller **SPOTSIZE**) and bring the image into **focus**).

- 2. Make sure the camera is **REMOVED** firstly and go to diffraction mode (**DIFF/IMG**). Insert and center the first **objective aperture (OA)**. The handling of the aperture knobs is described in figure 4.3.
- 3. Go to image mode (**DIFF/IMG**) and carefully insert the camera in «rabbit mode». Take an image with around 6000 counts. Compare the taken image with centered OA and the according image from part a) in focus. Describe the difference between them.
- 4. For your documentation: Explain these differences from what you have learned about the beam paths in a TEM in the course preparation. You can take an additional image with the smallest OA centered in diffraction mode.
- e) Magnification
- 1. This experiment can be included into experiment 1.3.
- 2. If not already done in the previous step, change the magnification of the microscope and take images of at least two different magnification settings. Always remove the camera, if you switch to lower magnification settings!
- 3. For your documentation: Describe what happens when you switch between different magnifications and explain your observation.

Magnification calibration with a cross grating (optional)

- 1. Follow this procedure if there is time left at day 1. Otherwise the instructor will give you images for the magnification calibration. You will need these to measure distances in your images and add scalebars to them.
- 2. Turn off the filament and let the instructor show you, how to change the sample holder.
- 3. Insert the sample holder containing the **cross-grating replica**.
- 4. Let the instructor check the alignment of the microscope. Take a look at it since you will do this from now on.
- 5. You should see a cross-grating (2160 lines/mm) with latex spheres ($\emptyset 0.261 \,\mu$ m).
- 6. Take images of the cross-grating at each (high mag) magnification setting (3 kx to 250 kx).

- Make sure the sample is in focus. Remember how you determined that the sample is in focus.
- At higher magnifications, make sure to get as many straight lines as possible on the camera (use the «rabbit mode» if necessary). Sometimes there are irregularities in the cross-grating that will make it hard to get a good calibration. Also try to get good images of the spheres to check your calibration.
- You will have to adjust the **spotsize** and **exposure time** as you go to higher magnifications.
- 7. Talk with the instructor how to measure the pixel size from the real space and the power spectrum (FFT) of the image using Fiji.
- 8. Understand how the power spectrum is calculated and qualitatively describe the differences in the power spectrum at lower and higher magnification.
- 9. You should now have at least one image of the cross-grating and latex spheres for each (possible) magnification.

Finishing day 1

- 1. Make sure you have recorded and saved all images for your documentation.
- 2. When done remove all apertures and close V3.
- 3. Remove the camera from the beam path.
- 4. Turn off the filament and then the high tension.
- 5. Remove the sample from the beam path.
- 6. Insert the sample you will investigate tomorrow (Au thin film).
- 7. When you have saved all your images (use a clean USB Stick), shut down the camera computer.
- 8. Turn off the microscope in the microscope software. (Make sure the camera computer has shut down completely.)
- 9. Shut down the microscope computer.
- 10. Finish Day 1.



Figure 4.2: Control panel for the ZEISS EM10EL.

4.2 Day 2: Diffraction Experiments

Firstly you will learn to align the microscope, since this is the everyday procedure in electron microscopy. Afterwards you will get to know electron diffraction and understand, how different apertures can be used for material investigation.

- 1. Turn on the microscope as described in the according section in the **«Alignment Procedure EM109EL»** document.
- 2. You can use the time the system needs for pumping to talk with the instructor about electron diffraction before you start over.
- 3. You will be measuring at 80 kV acceleration voltage and No. «2» emission settings in the software (around 5 to $6 \mu \text{A}$ emission current) unless otherwise stated.
- 4. Make sure no apertures are inserted.
- 5. Use the **«Alignment Procedure»** to align the microscope and record background and flatfield images for the camera.
- 6. Insert the first condensor aperture as described in the **«Alignment Proce**dure».

Get to know diffraction

- 1. Make sure the Au sample is inserted to the microscope.
- 2. Enter a small magnification in low-mag-mode and screen the sample to find an area containing the sample film.



Figure 4.3: (Condensor) aperture drive. Turning 1 will let you insert or remove the different sized apertures. The black knobs (2 and 3) let you move the aperture in x and y direction.

- 3. Increase the magnification of 20 kx.
- 4. Now make sure, that the camera is **NOT** in the beam path before you change to diffraction mode. You should look at the image on the fluorescent screen first.
- 5. Go to diffraction mode (DIFF/IMG).
- 6. Does the observation meet your expectation?
- 7. Go back to image mode (**DIFF/IMG**).
- 8. Insert and center the smallest selected-area aperture (SA).
- 9. Go back to diffraction mode (**DIFF/IMG**).
- 10. Describe and explain the differences.
- 11. Insert the biggest **objective aperture** (OA) and cover the central beam, so that you can use the camera in «rabbit mode».
- 12. Activate the «rabbit mode» and carefully insert the camera into the beam path. Make sure the central beam is still covered!
- 13. While in «rabbit mode» **move** the inserted **SA** back and forth in one direction while you are in diffraction mode and observe the diffaction pattern (DP) on the camera. **Describe** significant differences in the DP between two different SA positions.



Figure 4.4: (A) Air lock. Top Arrow indicates direction to turn when opening the air lock. Bottom arrow indicates the slider which is in parking position on this picture. (B) Specimen holder and tools. (C) Specimen holder when unscrewing the cap.

- 14. **Take** at least two images of the visible DP part at two different SA positions. It is <u>not</u> neccessary the reach 6000 counts to record a DP. Just check for a godd signal to noise ratio.
- 15. If done, remove the camera and switch to another SA of bigger size and center it.
- 16. As before **take** an image of the visible DP part with a bigger centered SA.
- 17. Explain your observations. Argue, what the selected-area aperture (SA) is useful for if you consider to record DPs.

Experiment 2.1: Investigating the diffraction pattern of a thin gold film

Record a full diffraction pattern (DP) using the **largest OA** at a camera length (CL) of 150 mm or 250 mm.

1. First decide, if it is reasonable to use a **SA** before recording the DP.

- 2. IMPORTANT!!! Make sure the (000) spot of the DP is NOT on the camera, but hidden by the OA.
- 3. IMPORTANT!!! Always remove the camera and use the fluorescent screen when adjusting the OA.
- 4. Use a reasonable **SPOTSIZE** and make sure your DP is in focus.
- 5. Move the **OA** around the center of the diffraction pattern to record a full DP of Au at the chosen camera length (CL). You will need about 6 to 8 images.
- 6. In your analysis you will load them as an "image sequence" into Fiji to prepare it for your documentation later on.
- 7. Remove the camera if you do not need it any more.

For the next day you have to decide now, which experiments you want to do since this will influence if you need to change the sample at this stage or not.

SEM: If you choose this option, proceed with experiment 2.2 and afterwards experiment 2.4. On the next day you will observe biological samples in the SEM.

TEM: If you choose this option, proceed with experiment 2.3 before experiment 2.2, so that you don't need to insert the Au film again. If there is no time left for experiment 2.3 today, you can perform it tomorrow. Skip experiment 2.4.

In the following experiment you will learn, how the OA can be used for darkfield imaging.

Experiment 2.2: Bright- and darkfield imaging

- 1. Remove the camera and turn off the filament to insert the MgO sample.
- 2. Use the **«Alignment Procedure»** to make sure the beam is aligned.
- 3. As in previous experiments investigate the sample in low-mag mode. **De-**scribe, what you see.
- 4. Toggle magnification in high-mag-mode, focus and spotsize to investigate your sample using the camera in «rabbit mode».
- 5. Find a sample area containing just a few MgO crystals, if possible.
- 6. Go to 20 kx or 30 kx and make sure your sample is in focus.

- 7. Insert and center the smallest SA on the camera.
- 8. When done, remove the camera and switch to diffraction mode (**DIFF/IMG**). You should see the DP on the fluorescent screen.
- 9. Insert and center the smallest **OA** (2nd aperture position).
- 10. Switch back to image mode (**DIFF/IMG**) and acquire a bright-field (BF) image. Make sure it is in focus (use the «rabbit mode» if necessary).
- 11. In diffraction mode center the OA around a diffracted beam. If you cover the central beam with the OA you can use the camera in «rabbit mode» to do fine adjustments. Take an image of the diffracted beam you used for your documentation.
- 12. Go to image mode (**DIFF**/**IMG**) and acquire a so called dark-field image.
- 13. Do this for a reasonable number of diffraction spots, and if possible of two opposing diffraction spots (Friedel pair). Feel free to go to higher magnifications or another sample position.
- 14. You should have acquired at least one BF image of a sample area and a few DF images from opposing and non-opposing diffraction spots. Later in your documentation you should be able to explain and discuss your observations.
- 15. Remove the SA and the OA.

Elective experiments

Choose one of the following experiments:

Experiment 2.3: Record another DP of the thin Au film at 150 mm or 250 mm camera length depending on which camera length you chose previously. You can proceed as described in experiment 2.1. This you will need later on to calibrate both camera lengths of the TEM.

Experiment 2.4: Choose at least one of the provided biological samples and insert it to the TEM. Take images of the regions of interest (e.g. amoeba and virus). On the next day you will compare them with images taken in the SEM.

Finishing day 2

- 1. When you finished all experiments and saved all images for your documentation, remove all apertures and close V3. Turn off the filament and then the high tension.
- 2. Remove the sample and the camera from the beam path.
- 3. Shut down the camera computer.
- 4. If neccessary (by choosing experiment 2.4) insert a sample you need on the next day. If you didn't have the time for experiment 2.3 enter the MgO sample otherwise choose either Au/Pd or Pd/Pt. Note which one you chose.
- 5. Turn off the microscope in the microscope software. (Make sure the camera computer has shut down completely.)
- 6. Shut down the microscope computer and turn off the main switch afterwards.
- 7. Finish day 2.

4.3 Day 3: Elective Experiments at the TEM or SEM

4.3.1 Elective Experiments at the TEM

- 1. Turn on the Microscope.
- 2. Use the **«Alignment Procedure»** to align the microscope and record background and flat-field images for the camera.

Experiment 3.1: Investigating Vegard's Law

- 1. Make sure, you inserted either a Au/Pd or Pd/Pt thin film into the microscope.
- 2. Record a full diffraction pattern (DP) as you did for the Au gold thin film in experiment 2.1 at camera lengths 150 mm and 250 mm.
- 3. Do you see a significant difference between the DPs of Au and the sample you chose?
- 4. When done, you should have recorded a set of images to create a full DP for either Au/Pd or Pt/Pd at camera lengths 150 mm and 250 mm for your documentation.

Experiment 3.2: Investigating Catalase Crystals

Similar to the cross-grating replica catalase is a standard for magnification calibration. You can use it to calibrate the high magnifications of the TEM, because the lattice spacings of catalase are much smaller. More interestingly it is possible to see the effect of the contrast transfer function at different defocus much better compared to the investigation of the myelin sheath of a mouse muscle.

- 1. Remove the camera and turn off the filament to insert the sample holder containing the **catalase sample**.
- 2. While you change the sample, explain the purpose of negative stain.
- 3. Once the filament has ramped up, change the emission setting to «3» in the software. You should now have a much larger emission current.
- 4. Use the **«Alignment Procedure»** to make sure the beam is aligned. After changing the emission current **new background and flat-field images** are needed for the camera.
- 5. Investigate the sample and describe what you see. You probably have to increase the magnification to see the crystal spacings of catalase. Find a catalase crystal.
- 6. Use the **OA** to increase the contrast.
- 7. Use the support film to correct the **objective astigmatism** (**Obj Stig**) at high magnification for best results. You might need to use a higher spotsize, longer exposure times and a sufficient high defocus to see the «Thon Rings» in the power spectrum. If the «Thon Rings» appear round, the objective astigmatism is minimized.
- 8. Choose a (high) magnification where you can see the «Thon Rings» and the crystal spacings in the power spectrum as well.
- 9. Change the focus to higher defocus and watch the crystal spacings and the power sectrum. **Describe** what you see.
- 10. Watch the crystal spacings until you realize changes in the image.
- 11. Observe the power spectrum closely and describe what happens there as well while changing the focus.
- 12. Record a **defocus series** as you already did in experiment 1.1. You may have to use another focus range!

- Start at focus and go to higher underfocus (defocus) in order to catch the changes in the images.
- Take images of your observations.
- 13. Later on in your documentation you should be able to discuss your observations.

Magnification calibration with catalase

- 1. Use the opportunity to take images of the catalase crystal to calibrate the higher magnifications of the microscope.
- 2. Bring the catalase crystal into focus. (use the «rabbit mode» if necessary).
- 3. Start at a magnification where you first see the catalase crystal spacings (note the power spectrum). Record images of the catalase crystal spacings. Do this for all higher magnification settings.
- 4. When done, you should have defocus series of the catalase crystal and images of it for each of the magnification settings in which you can see the spacings.
- 5. Use these images later in your documentation to calibrate the higher magnification settings. Using catalase the calibration can only be done from the Power Spectrum of the images.

Finishing Day 3

- 1. Make sure you have saved all images you need for your documentation and entered the microscope parameters in the excel sheet.
- 2. Go back to emission setting No. «2» (5 to $6 \mu A$), turn off the filament and afterwards the high tension.
- 3. Remove all apertures and close V3.
- 4. Remove the sample holder.
- 5. When you have saved all your images, shut down the camera computer.
- 6. Turn off the microscope in the microscope software. (Make sure the camera computer has shut down completely.)
- 7. Shut down the microscope computer.

- 8. Make sure you ask the instructor, if someting was unclear.
- 9. Finish Day 3.

4.3.2 Elective Experiments at the SEM

The Scanning electron microscope (SEM) is located in the BioQuant building 267.

- 1. The microscope will be turned on and adjusted by the instructor in the beginning. Let the instructor show you, how to toggle focus, magnification, stigmators and the apertures.
- 2. Get used to the controls and start optimising an image at a programmed sample position. You can always ask the instructor for help.

Experiment 3.4: Comparison between TEM and SEM

- 1. Move the sample holder to the samples that you already investigated in the TEM. You may have to search a bit for the regions of interest.
- 2. Since you don't want to damage the region of interest, choose a position next to it, to optimize the image settings.
- 3. Iteratively wobble the image and correct for astigmatism to get best results.
- 4. When done, take a reasonable amount of images and save them. You don't need to enter the microscope parameters in an excel sheet since they are saved in the images themselves.

Experiment 3.5: Investigating biological samples in the SEM

- 1. Move the sample holder to the sample position of other biological samples and take nice images.
- 2. Describe some of the problems you are facing while image acquisition and explain, why they occur. You should discuss this in your documentation.

5. Expected Documentation of the Experiments

Important! All real space images should contain a scalebar. In order to do this you need to calculate the pixelsize for the specific image firstly. Afterwards you can add a scalebar using Fiji (appendix B).

Indicate on your diffraction patterns the used camera length.

5.1 Magnification Calibration

- Create a calibration table containing the following for all magnification settings.
 - nominal magnification
 - pixel size and its error estimation
 - real magnification and the resulting error
 - method used for the calibration (real or fourier space)
 - the sample you used (cross grating or catalase)
 - check using the latex beads
- The neccessary information about the standard samples can be found in Appendix C. The course preparation from the beginning may help you here.
- Appendix B.3 explains how to calculate the power spectrum in Fiji.
- Plot the real magnification from the calibrated pixel size(detector pixel size: $14 \,\mu$ m) for each calibration method along with the nominal magnification. Compare real and nominal magnification of the microscope by a suitable fit function.
- Discuss the results and possible errors. Argue, why you need different specimens for calibrating all magnification settings by showing suitible images of the standard samples.
- Use the calibration to print a scalebar in all of the following images and to measure distances in real space and the power spectrum. Appendix B.1 explains, how to do this with Fiji. Make sure the scalebars are big enough.

5.2 Judging Image Quality

- Give a short explanation about the formation of Fresnel Fringes in your images. (No formulas!)
- Explain the effect of over- and underfocus using images of a holey carbon film. Describe, how the contrast of an image in focus can be increased. Figure 3.1 from your preparation may help you, since this figure was discussed in detail during the whole course.
- **Demonstrate the effect of astigmatism** on an image you recorded of a holey carbon film. A small schematic drawing may help here as well.
- Show, how coherence affects the image quality by judging the visibility of Fresnel diffraction patterns inside a hole in a carbon film.
- Give a short explanation, why the image rotates if you switch between the magnifications.

5.3 Diffraction, Bright- and Darkfield Imaging

- Explain diffraction shortly. Take the Bragg and Laue condition into account and show, how they are related.
- Explain the effect of an inserted selective aperture on the diffraction pattern (DP). You should have at least three images showing significant differences.
- Load the images from the DP as a stack into Fiji and use the Z-project function, to create a nice image of a full DP. How this can be done is explained in appendix B.1 and B.4.
- Index the diffraction pattern of gold. No calculation is needed here! (See the table A.1 in the appendix.)
- Explain the contrast differences between bright- and dark-field images from the MgO crystals. (especially the Friedel pair)
- Mention and discuss all other anomalies you may see in the DF images.

5.4 Understanding the Contrast Transfer Function

- Measure the spacings between the membranes in the myelin sheath using the power spectrum of the images. Appendix B.3 explains how to calculate the power spectrum of an image in Fiji.
- Using the power spectrum discuss, whether the measurement is correct. Take images from the SEM measurements into account, if you did experiment 3.4.
- If you consult the literature, you would find out, that there is no fix reference value for the distances in the myelin sheath. Discuss, why you would actually expect this.
- If you performed the elective experiment 3.2, show images of a Catalase crystal prepared as negative stain sample, and explain the purpose of «negative stain».
- Show the «Thon Rings» in the power spectrum of the myelin sample (and optional the catalase image) in over- and underfocus.
- Show and explain the influence of the contrast transfer function using the defocus series of the myelin in a mouse muscle (and optional the catalase crystal). (No formulas, but you can use a digram describing the CTF, if you find one!)

5.5 Investigating Vegard's Law

This section is only relevant, if you did experiment 3.1.

- Create a nice image of the diffraction pattern of Pt/Pd or Au/Pd using Fiji. (Appendix B.4)
- Calibrate the camera lengths at 150 mm and 250 mm using the according diffraction pattern of gold and Bragg's law. You can use the radial profile in Fiji for the calibration measurement (B.5).
- Analyze the diffraction patterns of Pt/Pd or Au/Pd at the camera lengths 150 mm and 250 mm respectively. From the rings in the DP you can measure the crystal spacings using Bragg's law and the calibration of the camera length. Don't forget to indicate the errors rising from your measurements.
- Calculate the lattice parameter for the sample you chose using Vegard's law and compare them with your measurements (Use Bragg's law). The ratio used to grow the cystals can be found in the appendix C.
- Discuss possible errors.

5.6 Comparison between TEM and SEM

This section is only relevant, if you did experiment 3.4 and 3.5.

- **Discuss** the advantages and disadvantages of imaging with TEM and SEM respectively. Discuss especially, why some samples can not be investigated in the TEM, but in the SEM.
- Show and discuss the images from your SEM measurements especially in comparison to your TEM investigation.
- Explain the occurence of sample charging in electron microscopy and how this influences your image.
- Discuss, why you didn't observe charging in the TEM.
- Discuss how you can avoid charging.

Bibliography

- [Ern10] R. ERNI. Aberration-corrected imaging in transmission electron microscopy: an introduction (ICP/Imperial College Press; Distributed by World Scientic Pub. Co, London : Singapore; Hackensack). 2010.
- [Hai98] M. HAIDER, H. ROSE, S. UHLEMANN, E. SCHWAN, B. KABIUS und K. URBAN. Aspherical-aberration-corrected 200 kV transmission electron microscope. Ultramicroscopy, 75, 1, 53–60. 1998.
- [Rei93] L. REIMER. Image Formation in Low-Voltage Scanning Electron Microscopy (SPIE - The International Society for Optical Engineering). 1993.
- [Rei08] L. REIMER und H. KOHL. Transmission Electron Microscopy, volume 36 of Springer Series in Optical Sciences (Springer New York). 2008.
- [Rod18] J. RODENBURG. The beginner's guide to transmission electron microscopy. URL http://www.rodenburg.org/guide/index.html. 2018.
- [Wil09] D. B. WILLIAMS und C. B. CARTER. Transmission Electron Microscopy: A Textbook for Materials Science (Springer Science & Business Media). 2009.
- [Wul16] S. F. WULF, V. ROPARS, S. FUJITA-BECKER, M. OSTER, G. HOFHAUS, L. G. TRABUCO, O. PYLYPENKO, H. L. SWEENEY, A. M. HOUDUSSE und R. R. SCHRÖDER. *Force-producing ADP state of myosin bound to actin.* Proceedings of the National Academy of Sciences, **113**(13), 1844– 1852. 2016.

Bibliography

Appendices

A. Diffraction Gold

h	k	1	$\mathrm{d}^* \; [\mathrm{nm}^{-1}]$	d [nm]
0	0	0	0	
1	1	1	4.256	0.235
0	0	2	4.914	0.204
0	2	2	6.949	0.144
1	1	3	8.149	0.123
2	2	2	8.511	0.117
0	0	4	9.828	0.102
1	3	3	10.71	0.093
0	2	4	10.988	0.091
2	2	4	12.037	0.083

 Table A.1: Peak positions for gold (Au).

A. Diffraction Gold

B. Helpful Features in Fiji

Fiji (Fiji is just ImageJ) is an opensource software for scientific image processing. You can use it to calibrate and analyze electron microscopy images. Fiji (including the Radial Profile Plugin) is installed on the CIP-Pool computers. Download sources and further information can be found on the internet:

• Fiji can be downloaded here for all common used operating systems:

http://imagej.net/Fiji/Downloads

• The Radial Profile Plugin here:

https://imagej.nih.gov/ij/plugins/radial-prole.html

• An extensive documentation for ImageJ in general can be found under:

https://imagej.nih.gov/ij/docs/guide/index.html

B.1 Opening images and scalebars

Single images can be open by drag and drop. You can open several images as one stack under:

File \rightarrow Import \rightarrow Image Sequence.

Fiji cannot read any calibration done by this camera software. Reset the image size to pixels for each image under:

Analyze \rightarrow Set Scale \rightarrow button «click to remove scale»

Here you can enter the results from your calibration as well, by entering the calculated pixel size in nm, once you know it. A scalebar can be added to your image under:

Analyze \rightarrow Tools \rightarrow Scale Bar

B.2 Adjusting Brightness and Contrast

The images you acquired during the experiment will be **16-bit greyscale**. Sometimes, especially at higher magnications, the intensity during acquisition was low. This might cause your image to appear black at first, since Fiji will use the whole (gray scale) histogram to display the image. You can adjust the histogram under:

Image \rightarrow Adjust \rightarrow Brightness/Contrast

Using the shortcut (Ctrl + C) works even quicker. Pressing the (Auto)-button in the window often yields to a nice view. After you have adjusted this to your liking, you can change the type of your image to 8-bit (Image -Type -8-bit). Fiji will then discard the information that you did not select in the histogram. Changing your images to 8-bit is sometimes helpful, since they will then take up less disk space.

B.3 Fast Fourier Transforms (Power Spectrum)

If you need to calculate the power spectrum of an image (e.g. for calibration), this can be done under:

 $\text{Process} \to \text{FFT} \to \text{FFT}$

This works even for a selected area of your image. When hovering the mouse pointer over a pixel of interest in the FFT, Fiji will display its position in the FFT and the respective distance in the real image. This can help you calibrate the magnification.

B.4 Image Stacks

Image stacks can be created from sequential images, while the slider in the window of an image stack represents the «Z-axis» of the stack. You can create a projection of all images in the stack by using:

Image \rightarrow Stacks \rightarrow Z-Project

When putting together the recorded partial diffraction patterns, choosing «Max Intensity» will let you create a single image containing the brightest parts of each single image. This is useful for creating a nice full diffraction pattern from your partial ones.

B.5 Radial and Line Profile

For your calibration you need to measure lengths in your images (e.g. cross-grating). For proper measurements (including error estimation) you should use «Line Profiles». Draw a line across the image using the «line tool». Double clicking the tool button lets you change the width of the line. You get a line profile under:

Analyze \rightarrow Plot Profile

Clicking the «List» button in the profile window will give you the x and y coordinates. You can export them (e.g. as .txt-file) and plot/analyze the profile in a software of your choice. The «live» button in the profile window allows you rearrange the drawn line in the image, while keeping the plot profile updated in real time. This is especially helpful for the radial profile.

The radial profile is similar to the plot profile. It measures the mean intensity of concentric circles as a function of the radius. Once you have installed the «Radial Profile» plugin, fit a circle to the data (e.g. of diffraction patterns). Make sure you press «Shift» so your circle becomes round and not elliptic. It is important that the plugin knows where the center of the circle is located since this will define the zero position of the radial profile. Be precise with fitting the circle. Otherwise the error of the measurement increases. Make use of the «live» button to optimize the circle position.

C. Provided Samples

- Inorganic samples
 - Gold (Au)
 - Magnesium oxide crystals
 - Gold/Palladium (Au/Pd 80:20)
 - Platinum/Palladium (Pt/Pd 80:20)
- Biological samples
 - Mouse muscle
 - Infected amoeba
 - Wasp (SEM)
 - Butterfly (SEM)
 - Pollen (SEM)
- Standard samples
 - Cross-grating calibration sample with latex spheres (2160 lines/mm, \varnothing 0.261 $\mu \rm{m})$
 - Bovine catalase